## Gebrauchsanleitung TPP Clipmax

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## Instructions for Use TPP Clipmax

*Clipmax is a combination of a cell culture flask and a slide made of cycloolefin polymer (COP) with high optical quality. Adherent cells can be cultivated on a growth area of 10 cm<sup>2</sup> and directly fixed and stained for microscopic examination. COP has excellent optical properties that allow measurements < 300 nm with a very low auto fluorescence. The good chemical resistance allows staining techniques with organic solvents.*

*For optimal cell adhesion and growth, only the slide is opto-mechanically activated.*

*The media chamber and slide are connected with a biocompatible seal. The chamber can be removed without tools by simply clipping it off. The system allows direct gas exchange through the filter screw cap.*

*Clipmax is only available with a filter screw cap.*

*Clipmax is for single use only.*

### Note:

*Follow the national regulations for handling biological materials and wear the appropriate protective clothing.*

*During the work process, be sure to observe the rules of aseptic technique.*

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### Instructions:

- *Inspect packaging and product for integrity. Use only materials in perfect condition.*
- *Check the expiry date (EXP) on the label and packaging. Use only products with a valid EXP.*
- *Open the package in a sterile environment and remove a complete system for use.*
- *Open the bioreactor and fill it with the medium and inoculum according to your laboratory routine.*
- *To obtain uniform cell growth, the following points must be observed:*
  - *Resuspend the cells well in the medium to avoid cell clumping.*
  - *Avoid foaming during resuspension and seeding to avoid holes in the cell layer.*
  - *Low volume of medium leads to meniscus formation and thus to uneven cell growth (cell clumps at the edge). Please refer to the technical data for the optimal filling volume.*
- *Close the filter screw cap with ventilation holes. Continuous gas exchange occurs through the integrated 0.22 µm hydrophobic membrane. If the membrane becomes wet, gas exchange is temporarily reduced.*
- *Cultivate the cells according to the usual laboratory routine.*
- *Aeration cams on the top of the chamber allow for even heat distribution in the incubator, even when multiple Clipmax are stacked.*
- *Fix and stain directly on the slide according to the protocol.*
- *Store at room temperature, protected from UV light.*

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**Remarks and Recommendations:**

- *Clipping the chamber from the slide is irreversible and results in a leaky chamber system.*
- *Use the labeling area on the Clipmax slide for identification.*
- *The stacking rim allows stable stacking of several Clipmax.*
- *For homogeneous cultivation of cells, it is essential to level the incubator and provide a continuous supply of medium.*
- *Avoid and/or eliminate sources of vibration near the incubator to ensure uniform cell growth.*

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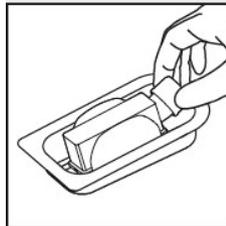
**Additional:**

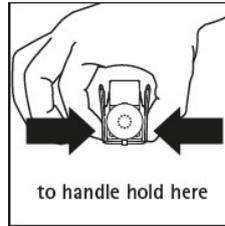
*Instructions for use, chemical resistance lists, and quality certificates for each product are available for download from [www.tpp.ch](http://www.tpp.ch).*

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