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### **TPP Dreifach-Verpackung 3-B SAL 10<sup>-6</sup>**

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## **Instruction for Use**

### **TPP Triple Bagged 3-B SAL 10<sup>-6</sup>**

*The 3-B packaging is specially developed for applications where sterile product transfer is essential. A high sterility assurance level (SAL) of 10<sup>-6</sup> ensures maximum sterility of the products.*

*3-B products are for single-use only.*

#### **Safety instructions:**

##### **1. Handling and Safety**

*When dealing with biological material, adherence to national regulations is mandatory. Furthermore, it is essential to ensure compliance with the established biological safety level of the laboratory, along with the safety data sheets and the manufacturer's instructions for use. Wear your personal protective equipment.*

##### **2. Risk of contamination**

*Always follow the rules for aseptic working and good laboratory practice. Inspect the packaging for integrity before use, as consumables are considered sterile only if the packaging is intact. Open the packaging immediately before use, and use only items that are visually flawless. Discard any products showing visible defects or signs of contamination.*

##### **3. Storage**

*Store TPP products at room temperature and protect them from direct UV radiation.*

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### **Instruction**

- *Check the expiration date (EXP) on the label and packaging. Only use products with a valid EXP date.*
- *Each product is triple-wrapped in three easy peel-open bags.*
- *For safety, the bag sizes are to prevent damage to the inner bag during opening, thus preventing contamination of the product.*
- *Each bag is marked with a security number:  
3 – outer bag,  
2 – middle bag,  
1 – inner bag.*
- *Bag 1 features product data: description, lot number, and EXP.*
- *Open the packaging in a sterile environment and remove the product.*
- *The product is now ready to be used for its intended purpose.*

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### **Usage and storage at sub-zero temperatures**

- **TPP Note:** Storage at subzero temperatures is not an intended application for this TPP product. TPP recommends transferring samples for these conditions into TPP cryo tubes, which are specifically designed for this purpose.
- Do not expose the consumables to liquid nitrogen. They may become damaged or burst when thawed.
- When using bioreactors or centrifuge tubes made of polypropylene (PP) at subzero temperatures, a preliminary application test should be conducted to ensure that the material withstands the specific requirements.

#### **Delivery time:**

Please note that this product may not be in stock but produced to order; a delay in delivery may occur.

#### **Additional:**

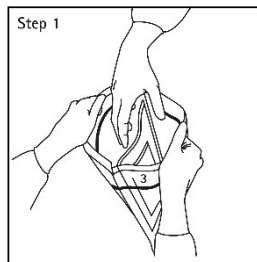
- SAL-value:  $10^{-6}$  for high sterilization requirements.
- Full TPP product line available worldwide.

Instructions for use, chemical resistance lists, and quality certificates for each product are available for download from [www.tpp.ch](http://www.tpp.ch).

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
**Step 1 / Bag 3,  
Entrance Area**

- Open bag 3 in the non-sterile environment.

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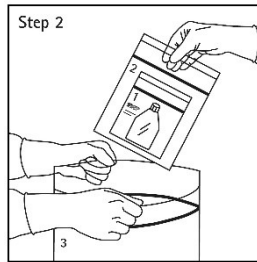
**Step 2 / Bag 2, Transfer**

- Pull out bag 2.
- Transfer to a cleanroom look.

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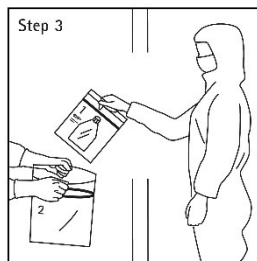**Step 3 / Bag 1, Cleanroom Look**

- Open bag 2 in the cleanroom look or sterile environment, and pull out bag 1.

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**Step 4 / Bag 1, Clean Room**

- Remove product from bag 1 under the biological safety cabinet.

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### **Disclaimer:**

*TPP products are intended solely for use in research and are not suitable for clinical, diagnostic, or therapeutic applications.*

*TPP assumes no liability for damages or defects resulting from improper or unauthorized use. Users are responsible for storing, handling, and using the products in accordance with the provided instructions.*

*TPP's recommendations are general guidelines and may not cover all possible scenarios. TPP is not liable for indirect, incidental, consequential, or special damages arising from the use or misuse of these products.*

*Swiss law governs these terms of use and any resulting legal matters.*