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Instruction for Use TPP Centrifuge Tube

The TPP centrifuge tubes are for preparing, mixing, centrifuging and storing suspensions (cells, bacteria, precipitates, etc.), solid and liquid reagents and samples.

The centrifuge tube is for single use only.

Safety instructions:

1. Handling and Safety

When dealing with biological material, adherence to national regulations is mandatory. Furthermore, it is essential to ensure compliance with the established biological safety level of the laboratory, along with the safety data sheets and the manufacturer's instructions for use.

Wear your personal protective equipment.

2. Risk of contamination

Always follow the rules for aseptic working and good laboratory practice. Inspect the packaging for integrity before use, as consumables are considered sterile only if the packaging is intact. Open the packaging immediately before use, and use only items that are visually flawless. Discard any products showing visible defects or signs of contamination.

3. Storage

Store TPP products at room temperature and protect them from direct UV radiation.

Instruction

- *Check the expiration date (EXP) on the label and packaging. Only use products with a valid EXP date.*
- *Before use, verify that the packaging is intact, as the consumable is only considered sterile if the packaging is undamaged.*
- *Open the packaging in a sterile environment and remove the product.*
- *Open the screw cap and fill the tube with the medium and inoculum according to the usual laboratory routine.*

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Notes and Recommendations:

- *Centrifuging at (too) high speed may cause deformation of the tube.*
- *The graduation serves as a reference for the fill volume and is not an absolute value.*
- *Alcohol (90%) combined with mechanical stress (rubbing) may dissolve or blur the white labeling.*

Usage and storage at sub-zero temperatures

- **TPP Note:** *Storage at subzero temperatures is not an intended application for this TPP product. TPP recommends transferring samples for these conditions into TPP cryo tubes, which are specifically designed for this purpose.*
- *Do not expose the consumables to liquid nitrogen. They may become damaged or burst when thawed.*
- *When using centrifuge tubes made of polypropylene (PP) at subzero temperatures, a preliminary application test should be conducted to ensure that the material withstands the specific requirements.*

Important:

- *Temperatures below 0 °C cause reduced elasticity of polypropylene. Do not subject frozen tubes to mechanical stress, such as dropping or hitting them.*
- *When storing at sub-zero temperatures, ensure that the contents freeze evenly and from bottom to top. → If the contents of the tube freeze first at the top, expansion cracks may occur at the bottom. Styrofoam racks have a strong insulating effect, which makes it difficult for the contents of the tubes to freeze evenly.*

Additional:

Instructions for use, chemical resistance lists, and quality certificates for each product are available for download from www.tpp.ch.

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Disclaimer:

TPP products are intended solely for use in research and are not suitable for clinical, diagnostic, or therapeutic applications.

TPP assumes no liability for damages or defects resulting from improper or unauthorized use. Users are responsible for storing, handling, and using the products in accordance with the provided instructions.

TPP's recommendations are general guidelines and may not cover all possible scenarios. TPP is not liable for indirect, incidental, consequential, or special damages arising from the use or misuse of these products.

Swiss law governs these terms of use and any resulting legal matters.