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## **Instructions for Use of TPP Vacuum Filtration «rapid»-Filtermax**

The TPP filtration system «rapid»-Filtermax is suitable for sterile filtration of cell culture media, sera and aqueous solutions.

The 0.22 µm polyether sulfone (PES) membrane ensures high flow rates with low protein binding and foam formation. The «rapid»-Filtermax is available as a filtration unit (bottle top filter incl. reservoir flask) as well as a separate bottle top filter or flask, with a DIN GL45 thread.

The filtration system is available in volume sizes 150, 250, 500 and 1000 mL. Each system component is identical in construction, available separately, and can be combined with each other.

The square foot of the reservoir ensures a safe stand during sterile filtration.

The «rapid»-Filtermax is for single use only.

### **Safety instructions:**

#### **1. Handling and Safety**

When dealing with biological material, adherence to national regulations is mandatory. Furthermore, it is essential to ensure compliance with the established biological safety level of the laboratory, along with the safety data sheets and the manufacturer's instructions for use.

Wear your personal protective equipment.

#### **2. Risk of contamination**

Always follow the rules for aseptic techniques and good laboratory practices. Open the packaging only immediately before use and use only items that are visually flawless. Dispose of any products that show visible defects or signs of contamination.

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### 3. Storage

Store the TPP products at room temperature and protect them from direct UV radiation.

#### Instruction

- Check the expiration date (EXP) on the label and packaging. Only use products with a valid EXP date.
- Before use, verify that the packaging is intact, as the consumable is only considered sterile if the packaging is undamaged.
- Open the packaging in a sterile environment and remove the product.
- Check the filtration system to ensure that the bottle top filter is securely attached to the bottle. Tighten it if necessary, but avoid overtightening to prevent damaging the thread.

#### Alternative:

- If you use only the TPP bottle top filter, screw it securely onto a suitable sterile bottle, taking care not to damage the thread.
- Connect the hose of the regulated vacuum source to the hose adapter. Then insert the adapter into the connection of the bottle top filter.
- Position and secure the filtration system to prevent it from tipping over during the process.
- Remove the lid from the bottle top filter and carefully pour in the solution to be filtered.
- Apply the vacuum and filter until the bottle top filter is empty or the reservoir / bottle is full.
- Turn off the vacuum after filtration, place the lid on the bottle top filter, and disconnect the hose.
- Unscrew the bottle top filter, unpack the individually wrapped sterile screw cap, and close the bottle with it.

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- *Before regular use of the product, perform a test run with the chosen settings to verify its suitability.*

#### **Advice and Recommendations:**

- *The volume scale is an indication of the filling volume, not an absolute value.*
- *Alcohol 90% with mechanical stress (rubbing) may dissolve or blur the white color of the print.*

#### **Information**

- *Provide the following materials: vacuum source; vacuum tubing.*
- *Centrifuge or prefilter solutions with high particle densities as a priority to prevent filter clogging.*
- *Attach solutions with a low particle density and a high filtration throughput to another reservoir for additional filtration. Please observe the rules of aseptic technique.*
- *Once the pouch is open, keep the product in a sterile area to ensure sterility.*
- *The flask is not suitable for freezing.*
- *Do not autoclave product.*
- *This product is not for direct patient care or diagnostic procedures, but for research use only.*

#### **To avoid possible injury due to implosions:**

- *Do not exceed 700 mm Hg differential vacuum at 25 °C*
- *Bottle top filter on alternative bottles: Use only sterile plastic or glass bottles that are approved for sterile filtration under vacuum. These should be free of any damage (risk of implosion). Observe the maximum total pressure.*

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- *Wear eye-protection and perform the filtration in special safety units.*
- *Do not use hand-held vessels while filtering under vacuum.*

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**Usage and storage at sub-zero temperatures**

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- **TPP Note:** Storage at subzero temperatures is not an intended application for this TPP product. TPP recommends transferring samples for these conditions into TPP cryo tubes, which are specifically designed for this purpose.
- *Do not expose the consumables to liquid nitrogen. They may become damaged or burst when thawed.*

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*Instructions for use, chemical resistance lists, and quality certificates for each product are available for download from [www.tpp.ch](http://www.tpp.ch).*

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Filter membrane Polyether sulfone (PES),  
 type TPP fast flow  
 Flask and bottle PS  
 top filter  
 Screw cap PE  
 Vacuum connector PP

| Technische Daten                  | Technical Data                     | Werte / Values                                  |
|-----------------------------------|------------------------------------|---|
| Max. Gebrauchs-Temperatur °C      | Max temp in use °C                 | max. 45   |
| Vakuum mm Hg bei 25 °C            | Vacuum mm Hg at 25 °C              | 700   |
| Totvolumen ml                     | Dead volume mL                     | 2.3   |
| Vakuumanschluss mm                | Vacuum connection mm               | 6 – 10  |
| Wasserflussrate bei 25 °C         | Water flow rate at 25 °C           | 22 ml/min/cm <sup>2</sup><br>0.07 mPa (0.7 bar) |
| Proteinbindung µg/cm <sup>2</sup> | Protein binding µg/cm <sup>2</sup> | < 20  |

|                              |                             | <b>System<br/>System</b> | <b>Flaschenaufsatzfilter<br/>Bottle Top Filter</b> | <b>Flasche<br/>Flask</b> |
|------------------------------|-----------------------------|--------------------------|--|--------------------------|
| <b>Produkt Nr.</b>           | <b>Product No.</b>          | <b>99150</b>             | <b>99155</b>                                       | <b>99157</b>             |
| <b>Abmessungen</b>           | <b>Measurements</b>         |                          |  |                          |
| Höhe mm                      | Height mm                   | 103                      | 57   | 63                       |
| Breite x Länge<br>mm         | Width x Length<br>mm        | 93 x 93                  | 90 x 89  | 93 x 93                  |
| Filtergrösse cm <sup>2</sup> | Filter size cm <sup>2</sup> | 49                       | 49   | ---                      |
| Porengrösse µm               | Pore size µm                | 0.22                     | 0.22   | ---                      |
| Volumen ml                   | Volume mL                   | 150                      | 150  | 150                      |

|                              |                             | <b>99250</b> | <b>99255</b> | <b>99257</b> |
|------------------------------|-----------------------------|--------------|--------------|--------------|
| <b>Produkt Nr.</b>           | <b>Product No.</b>          |              |              |              |
| <b>Abmessungen</b>           | <b>Measurements</b>         |              |              |              |
| Höhe mm                      | Height mm                   | 143          | 75           | 87           |
| Breite x Länge<br>mm         | Width x Length<br>mm        | 93 x 93      | 90 x 89      | 93 x 93      |
| Filtergrösse cm <sup>2</sup> | Filter size cm <sup>2</sup> | 49           | 49           | ----         |
| Porengrösse µm               | Pore size µm                | 0.22         | 0.22         | ---          |
| Volumen ml                   | Volume mL                   | 250          | 250          | 250          |

|                              |                             | <b>99500</b> | <b>99505</b> | <b>99507</b> |
|------------------------------|-----------------------------|--------------|--------------|--------------|
| <b>Produkt Nr.</b>           | <b>Product No.</b>          |              |              |              |
| <b>Abmessungen</b>           | <b>Measurements</b>         |              |              |              |
| Höhe mm                      | Height mm                   | 213          | 111          | 121          |
| Breite x Länge<br>mm         | Width x Length<br>mm        | 93 x 93      | 90 x 89      | 93 x 93      |
| Filtergrösse cm <sup>2</sup> | Filter size cm <sup>2</sup> | 49           | 49           | ---          |
| Porengrösse µm               | Pore size µm                | 0.22         | 0.22         | ---          |
| Volumen ml                   | Volume mL                   | 500          | 500          | 500          |

|                              |                             | <b>99950</b> | <b>99955</b> | <b>99957</b> |
|------------------------------|-----------------------------|--------------|--------------|--------------|
| <b>Produkt Nr.</b>           | <b>Product No.</b>          |              |              |              |
| <b>Abmessungen</b>           | <b>Measurements</b>         |              |              |              |
| Höhe mm                      | Height mm                   | 285          | 143          | 160          |
| Breite x Länge<br>mm         | Width x Length<br>mm        | 111 x 111    | 108 x 108    | 111 x 111    |
| Filtergrösse cm <sup>2</sup> | Filter size cm <sup>2</sup> | 69           | 69           | ---          |
| Porengrösse µm               | Pore size µm                | 0.22         | 0.22         | ---          |
| Volumen ml                   | Volume mL                   | 1000         | 1000         | 1000         |

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### **Disclaimer:**

*TPP products are intended solely for use in research and are not suitable for clinical, diagnostic, or therapeutic applications.*

*TPP assumes no liability for damages or defects resulting from improper or unauthorized use. Users are responsible for storing, handling, and using the products in accordance with the provided instructions.*

*TPP's recommendations are general guidelines and may not cover all possible scenarios. TPP is not liable for indirect, incidental, consequential, or special damages arising from the use or misuse of these products.*

*Swiss law governs these terms of use and any resulting legal matters.*