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Instruction for Use TPP Cell Scraper and Spatula (Lifter)

The TPP cell scraper and Spatula (or lifter) are used for harvesting adherent cell cultures in various containers such as cell culture dishes, plates and bottles.

The angled and pivoting blade simplifies the harvesting procedure and allows for complete cell harvesting with minimal cell damage.

The cell scraper and cell lifter is for single use only.

Safety instructions:

1. Handling and Safety

When dealing with biological material, adherence to national regulations is mandatory. Furthermore, it is essential to ensure compliance with the established biological safety level of the laboratory, along with the safety data sheets and the manufacturer's instructions for use.

Wear your personal protective equipment.

2. Risk of contamination

Always follow the rules for aseptic techniques and good laboratory practices. Open the packaging only immediately before use and use only items that are visually flawless. Dispose of any products that show visible defects or signs of contamination.

3. Storage

Store the TPP products at room temperature and protect them from direct UV radiation.

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Instruction:

- *Check the expiration date (EXP) on the label and packaging. Only use products with a valid EXP date.*
- *Before use, verify that the packaging is intact, as the consumable is only considered sterile if the packaging is undamaged.*
- *Open the packaging in a sterile environment and remove the product.*

Fields of Application:

- *The cell scrapers are available in 3 sizes to suit your needs (see selection guide).*
- *In addition to traditional cell culture applications, the sterile, individually packaged scrapers and spatulas (lifter) can be used for a variety of other applications.*

Additional:

Instructions for use, chemical resistance lists, and quality certificates for each product are available for download from www.tpp.ch.

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Disclaimer:

TPP products are intended solely for use in research and are not suitable for clinical, diagnostic, or therapeutic applications.

TPP assumes no liability for damages or defects resulting from improper or unauthorized use. Users are responsible for storing, handling, and using the products in accordance with the provided instructions.

TPP's recommendations are general guidelines and may not cover all possible scenarios. TPP is not liable for indirect, incidental, consequential, or special damages arising from the use or misuse of these products.

Swiss law governs these terms of use and any resulting legal matters.